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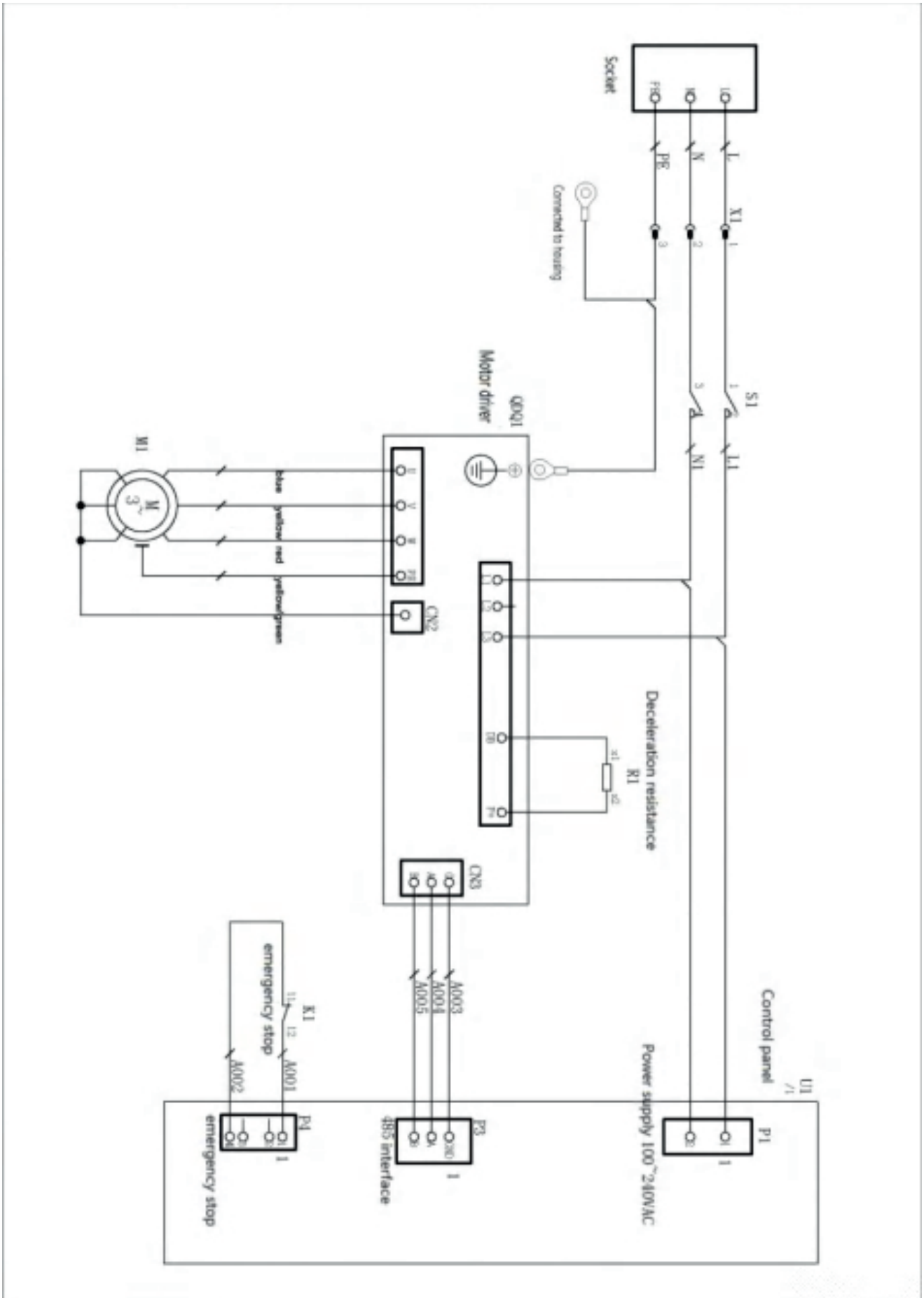
## SAFETY NOTICE

- » Please read this manual thoroughly before installing and using this product. Pay special attention to the content and recommendations, and follow the safety precautions to avoid equipment damage, fire, and personal injury.
- » Ensure that the grounding meets the required standards.
- » Operators must complete safety training before starting work; only qualified personnel may operate this machine.
- » Make sure the environment meets the installation requirements. Operate the equipment according to the operational guidelines.
- » Always disconnect the equipment from the power supply before cleaning.
- » In case of malfunction, do not disassemble the device. Instead, seek professional maintenance to avoid electrical shock.
- » Please keep this manual in good condition.

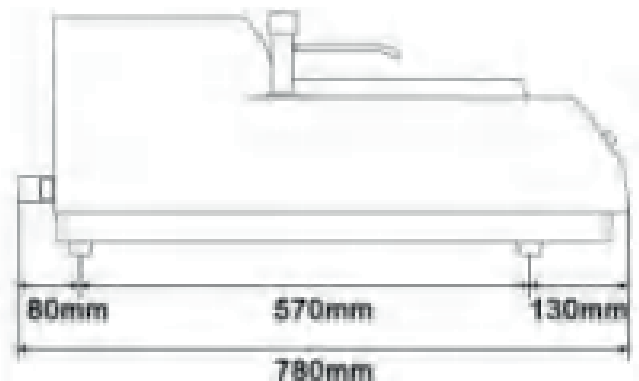
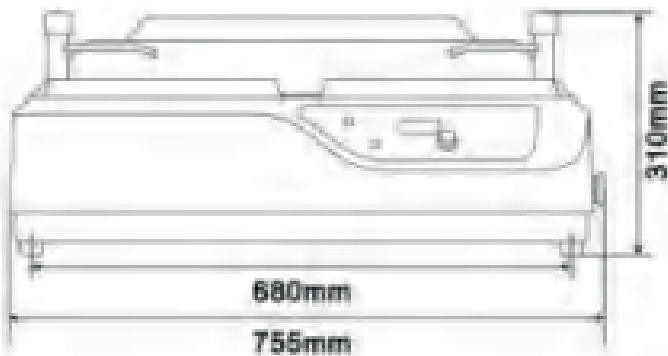
## INTRODUCTION

This grinding and polishing machine features two rotating wheels. It is suitable for pre-grinding, grinding, and polishing metallographic samples. The machine is controlled by a microprocessor, allowing for speed adjustments from 0 to 1400 RPM, ensuring a wide range of applications. The device is equipped with water cooling to cool the sample during grinding, preventing overheating and damage to the metallographic structure of the sample. With easy operation and reliable performance, this machine is an essential tool for factories, research institutions, and universities for preparing metallographic samples.


## TECHNICAL SPECIFICATIONS

- » Specification


<b>Disc Diameter</b>	Ø203 mm
<b>Disc Rotation Speed</b>	0 - 1400 RPM (stepless adjustable)
<b>Disc Rotation Direction</b>	Clockwise or counterclockwise
<b>Power Supply</b>	AC 220 V, 50 Hz
<b>Motor Power</b>	0.75 kW
<b>Dimensions</b>	755×700×310 mm
<b>Weight</b>	50 kg

## » Machine Diagram

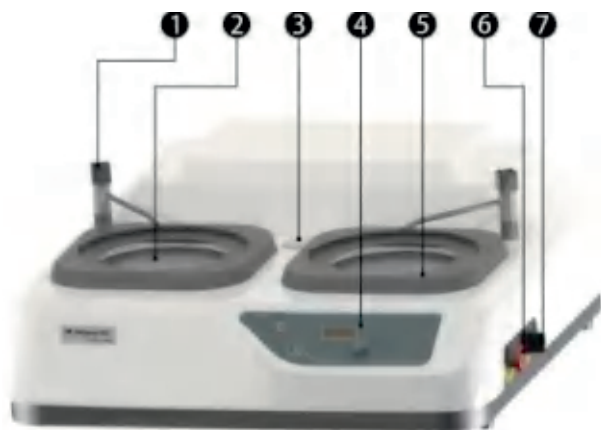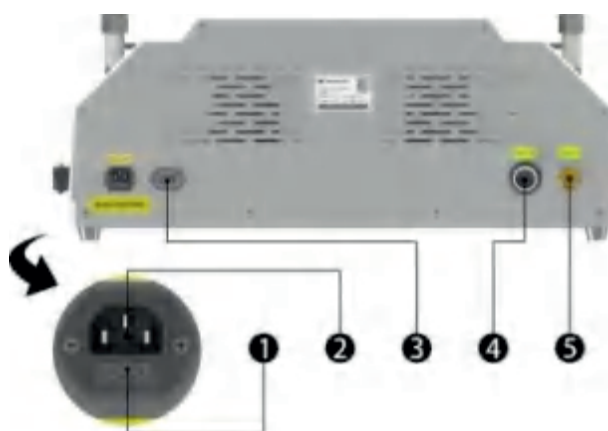- ① Water valve      ② Disc      ③ Head port
- ④ Control panel    ⑤ Locking ring
- ⑥ Emergency stop   ⑦ Switch

- ① Safety socket      ② Power cord connection
- ③ Head connection (head is optional)
- ④ Drain pipe connection    ⑤ Water inlet connection


## INSTALLATION AND PRECAUTIONS

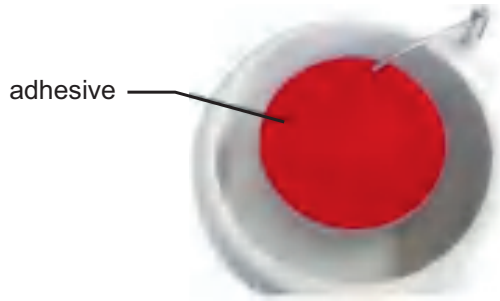1. The device must be properly grounded and should use a three-pronged power outlet that meets standard requirements.
2. The device should be installed on a stable workbench that does not produce vibrations. The environment should be ventilated, dry, with a temperature range of 10–30°C and relative humidity below 85%. There should be no corrosive gases or conductive dust.
3. Unpack the box, remove the accessories, and place the equipment on the workbench. Insert one end of the drain pipe into the outlet at the back of the device, ensuring it is securely fitted. Connect the water inlet to the corresponding input, making sure the rubber gasket contacts the inlet properly to form a tight seal.
  - ⚠ The drain pipe must be positioned lower than the device body to ensure smooth drainage without leaks.
4. Determine the device's placement based on the position of the water inlet and outlet. Ensure the device is level and stable.
5. Remove the locking ring from the device, take out the disc from the accessories, and clean the tapered area of the disc. While standing in the working position, clean the tapered surface of the main shaft inside the device, then install the disc onto the main shaft.
  - ⚠ Due to the connection between the disc and the main shaft, use a rubber mallet to gently tap the center of the grinding disc until it fits snugly. Improper installation may affect the machine's performance.

## 6. Metallographic Grinding Paper and Polishing Cloth Installation:

### 1) For PSA (Pressure-Sensitive Adhesive) Grinding Paper and Polishing Cloth:


Clean the top surface of the disc and attach the grinding paper or polishing cloth using special adhesive. Visually ensure the distance between the paper or cloth and the disc's edge is uniform. Press the paper or cloth evenly by hand to ensure firm adhesion.

### 2) For Non-Adhesive Grinding Paper and Polishing Cloth:

a) Clean the top surface of the disc, place the non-adhesive grinding paper or polishing cloth on the disc, and then secure it by placing the locking ring on the disc's outer edge.

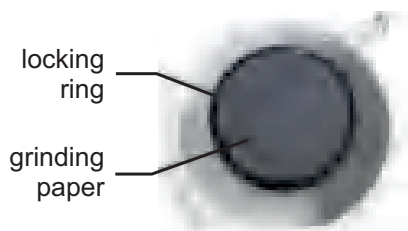
b) If using a microfiber polishing cloth, ensure it lies flat on the disc, with its edges evenly aligned. Then, place the locking ring on the disc's outer edge to secure the cloth.

7. After closing the switch on the external outlet, connect the water supply hose to the external outlet. Ensure the water valve is closed, then open it after connection. No water leakage should occur at the connection points.


8. Place the drain pipe into a container or connect it to a drainage system for waste liquid disposal. Open the water valve and check for smooth water flow through the inlet and drain pipes, ensuring no leaks at connection points. Then, close the water valve.

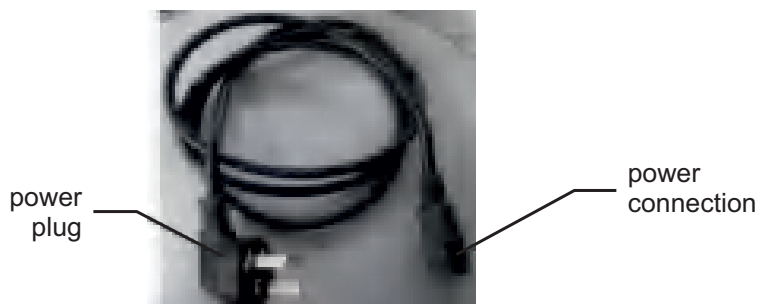9. Connect the power cord to the device.

## CONTROL PANEL AND DISPLAY

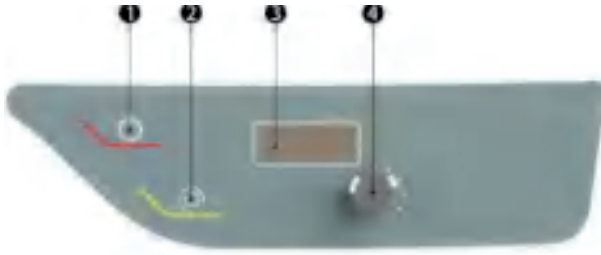The control panel is shown below:

1. Start/Stop: Press this button to start the disc; press it again to stop.
2. Control Key: Press to change the direction of the disc's rotation.
3. Display: Shows the set disc rotation speed.
4. Control Knob: Adjusts the disc rotation speed (RPM).

## DESCRIPTION OF OPERATION AND WARNINGS

1. Plug the power cord into the socket. The device will power on, and the display will show the set rotation speed. Press the start button to begin operation; the disc will start spinning. Allow the device to run without a load for approximately 1 minute to check for normal operation before proceeding to the next step.

⚠ Press the emergency stop button to halt the device if any danger arises during grinding or polishing.

⚠ The emergency stop button should not be used as a regular stop button.

2. Grinding Method:

After cutting, metallographic samples should be prepared using a grinder and cooled with water to avoid unwanted changes in the material's structure. Samples cut with a specialized cutting machine can be directly used for rough grinding.

Grinding is divided into three steps: rough grinding, semi-fine grinding, and fine grinding.

Table 1: Grinding parameters for the grinding and polishing block:

Type of Grinding	Grinding Paper Grit	Disc Speed	Requirements
<b>Rough Grinding</b>	240# - 400#	50 RPM - 1000 RPM	Remove machining marks from cutting.
<b>Semi-Fine Grinding</b>	600# - 800#	50 RPM - 1000 RPM	Remove marks from rough grinding.
<b>Fine Grinding</b>	1000# - 2000#	50 RPM - 1000 RPM	Achieve a near-mirror surface, removing visible marks.

Note: The abrasive size and disc speed should be selected based on the requirements of different metallographic samples. For specific samples, the semi-fine grinding step can be skipped. If a higher sample demand exists, additional grinding steps can be added as necessary.

### Grinding Method:

a) First, select an appropriate coarse metallographic grinding paper and install it on the disc. Press the start button to start the device and set the speed to 150 RPM; use the knob to adjust the disc to the desired speed. Slowly turn the control knob on the water valve to open it and set the water flow, adjusting from dripping to a steady stream as needed. Firmly hold the cut sample and gently press it against the grinding paper. Initially, apply light pressure in the center of the grinding disc while gradually moving the sample outward. Once the sample makes proper contact with the grinding paper without deviation, increase the pressure to approximately 2 kgf to prevent overheating and damage to the ground surface due to friction. Rotate the sample 90° with each change of grinding paper to thoroughly remove previous grinding marks until the requirements are met. Then press the stop button to end the operation and close the water valve.

△ Ensure proper drainage to prevent excessive water accumulation in the device, which could seep through the inner walls.

△ After pressing the stop button, wait briefly before pressing the start button again; the device will resume operation at the speed set before stopping.

b) For selected semi-fine or fine grinding, install the appropriate metallographic grinding paper on the disc. The grinding method is the same as for coarse grinding and continues until the fine grinding requirements are met. At the end of the work, press the stop button to halt operation and close the water valve, concluding the grinding process.

### 3. Polishing Methods:

Polishing consists of rough and fine polishing.

Type of Polishing	Abrasive Particle Size	Disc Speed	Requirements
<b>Rough Polishing</b>	W3 - W5	50 RPM - 1000 RPM	Remove fine grinding marks and surface layers.
<b>Fine Polishing</b>	W0.5 - W2	50 RPM - 1000 RPM	Achieve fine polishing and smooth, scratch-free surface.

#### Note:

1. The type of abrasive, particle size, disc speed, and polishing cloth should be selected based on the requirements of different metallographic samples. Common supplies for grinding/polishing include polishing pastes, sprays, and abrasive suspensions. Materials may include aluminum oxide, chromium oxide, diamond, and colloidal silica.
2. If the sample requirements are high, additional polishing steps may be added.

### Polishing Method:

a) **Rough Polishing:** Install adhesive or non-adhesive polishing cloth on the disc. Turn the control knob on the water valve to open it, allowing water to drip onto the cloth. Press the start button to begin operation. Set the device to run at 300 RPM; use the knob to adjust the disc to the desired speed. Once the cloth is fully moistened, close the water valve and apply the modified polishing agent to the cloth for approximately 10 seconds, pouring off excess water beforehand. Firmly hold the sample and gently press it against the cloth. Initially, apply light pressure at the center of the disc, gradually moving outward. Once the sample contacts the cloth without deviation, increase the pressure. However, excessive pressure should be avoided to prevent motor overload and cloth tearing. Rotate the sample 90° with each change of polishing cloth to remove previous marks thoroughly. If the cloth becomes very sticky during operation, dilute the polishing agent further until rough polishing requirements are met. Finally, press the stop button to halt operation.

- b) Finishing Polishing: Install adhesive or non-adhesive polishing cloth on the disc. The method for finishing polishing is the same as for rough polishing, using finer polishing abrasives until the sample meets metallographic analysis requirements. At the end of the work, press the stop button to halt operation; turn off the power switch, system power, and water supply.
4. Plastic components of this machine should be used at temperatures above 0°C with special care to prevent damage.
  5. If abnormal noises are detected from the device, stop the machine immediately for inspection.
  6. After each operation, clean and maintain the device.
  7. Regularly inspect the device's belt for looseness, aging, or cracks, as these issues can significantly affect its performance.
- ⚠ Do not use damaged grinding paper or polishing cloth, as this may pose a danger during grinding or polishing.

## TROUBLESHOOTING

Problem Type	Troubleshooting Methods
Display does not turn on after pressing the power button	<ol style="list-style-type: none"> <li>1. Check if the power is properly connected.</li> <li>2. Inspect the power cables and sockets for damage.</li> <li>3. Verify that the fuse is not blown.</li> </ol>
Disc does not rotate after pressing the start button	<ol style="list-style-type: none"> <li>1. Check if the control buttons are damaged.</li> <li>2. Inspect the motor for any faults.</li> </ol>
Noise of disc collision with piston shaft during startup	Replace the worn locking ring if it is damaged.
Insufficient performance after extended use	Check the tension and wear of the drive belt.

#	Displayed Abbreviation on Screen
1	OC1: Overload 1
2	OC2: Overload 2
3	EH: Sampling Circuit Damage
4	PLD: CPLD Error
5	OS: Overspeed
6	HU: Overvoltage
7	EC: Abnormal Communication with Encoder
8	EC2: Abnormal Communication with Encoder 2
9	EP: Production Circuit Error
10	AH: Drive Overheating

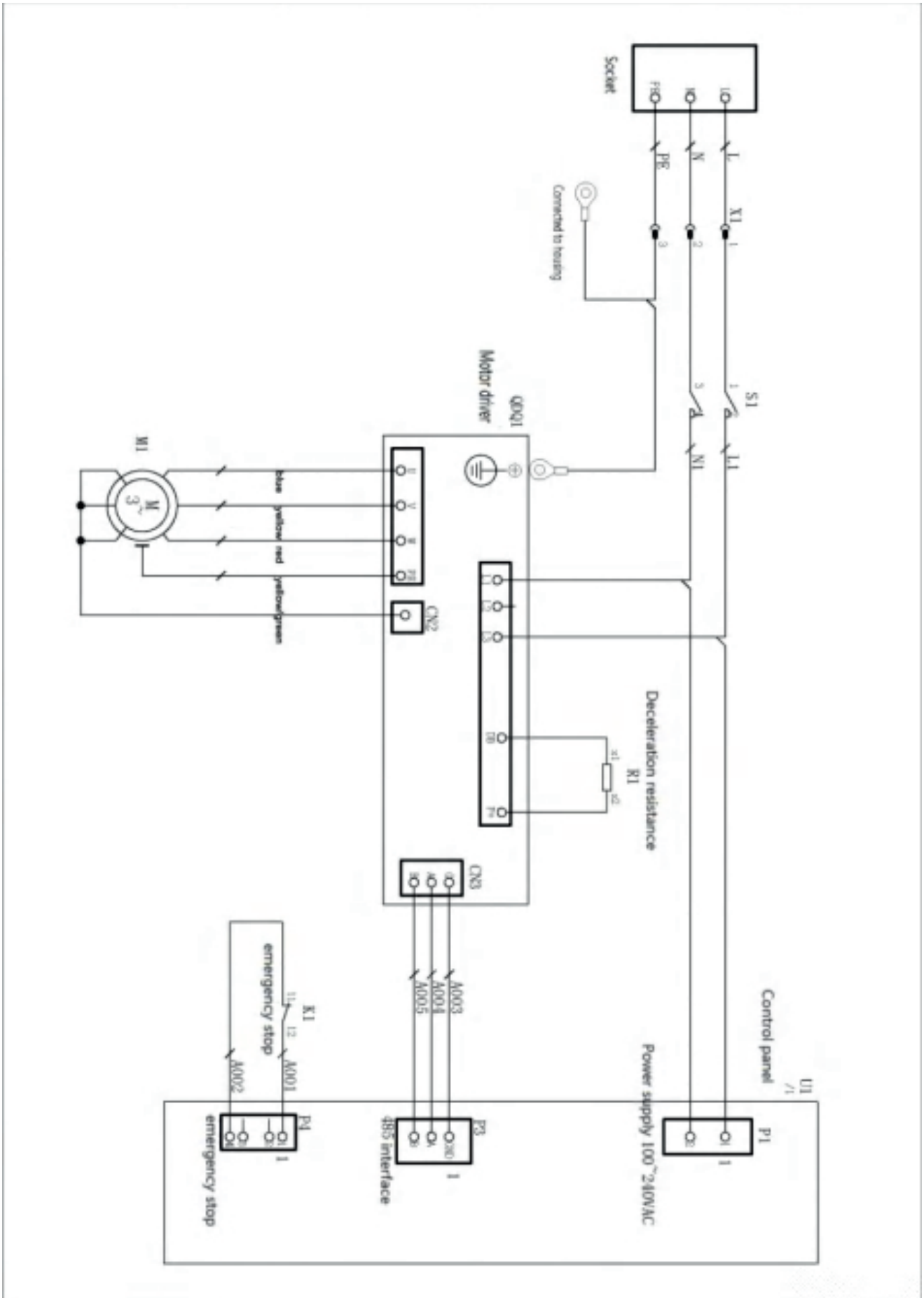
11	RH: Overheating of Regenerative Resistance
12	DE: Memory Exception
13	CE: Motor Code Error
14	SE: Overspeed Error
15	OL: Overload
16	OL2: Overload 2
17	LU: Low Voltage
18	OF: Deviation
19	POL: Power Outage Alarm
20	ND: Missing Motor Code
21	CE: Motor Selection Error
22	GOH: Zero Error
23	PLOT: Soft Limit Alarm
24	PNOT: Negative Soft Limit Alarm

Note: After resolving the issues, reset the device by turning off the power.

#	Displayed Abbreviation on Screen
1	E100: Stop Press
2	E101: Start Button Long Press
3	E102: Control Key Long Press
4	E103: Left Button Long Press
5	E104: Right Button Long Press
6	E105: V1 Button Long Press
7	E106: V2 Button Long Press
8	E107: V3 Button Long Press
9	E108: V4 Button Long Press

Note: After resolving the issues, press the emergency stop button and then release it to reset.

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**INSIZE is a global manufacturer of measuring technology** bringing technological innovations with representation in 75 countries around the world. Measuring instruments of the INSIZE brand represent the optimal solution without compromises to meet even the most demanding measuring needs.

**INSIZE measuring instruments will pleasantly surprise you with:**  
| high-quality craftsmanship | reliable performance | affordable prices

INSIZE provides a complete portfolio of measuring instruments » calipers, height gauges, dial indicators, micrometers, roughness testers, hardness testers, measuring microscopes, optical measuring devices, profile projectors, testing machines, contour gauges, roundness measuring machines, thickness gauges, gages, protractors, force gauges, meters, scales, videoscopes, torque wrenches and accessories including stands for measuring instruments.

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